

FEATURES

- High Performance Scientific Camera
 - Hamamatsu S10141 Sensors
 - 1024 x 122 to 2048 x 506 active pixels
 - 12 μm x 12 μm pixel size
- Integrated TE Cooling
 - -15 $^{\circ}\text{C}$ from 20 $^{\circ}\text{C}$ ambient
 - Hermetically sealed
 - Temperature controlled to ± 0.1 $^{\circ}\text{C}$
 - Air cooled
- Integrated Digital Signal Processing
 - Utilizes [MityDSP](#) technology
 - Can be customized for application specific processing
- Interface Options
 - USB 2.0
 - 100 Mbit Ethernet
- Programmable General Purpose I/O
 - Use as trigger inputs or drive as shutter / frame strobes

APPLICATIONS

- Embedded Instrumentation
- Low Light Imaging Applications
- Raman Spectroscopy
- Portable Scientific Instrumentation
- Astronomy



Figure 1: MityCCD-H10141 with Standard Body

DESCRIPTION

The MityCCD-H10141 family of cameras integrates thermoelectrically cooled high performance CCD sensors suitable for scientific instrumentation with digital signal processing (DSP) capabilities.

An embedded camera rich in features, MityCCD is ready for stand-alone applications. A simple block diagram of the MityCCD camera is illustrated in Figure 2.

By integrating a floating point DSP and field programmable gate array into the camera electronics, the MityCCD camera provides the capability to embed application specific code (binning patterns, post processing algorithms, etc.) into a delivered instrument.

The MityCCD uses standard high speed interfaces including 100 Mbit Ethernet and USB 2.0 and eliminates the need for custom interface cards or electronics to a display PC, allowing the use of laptop devices, embedded computers, or (via Ethernet) remote display machines.

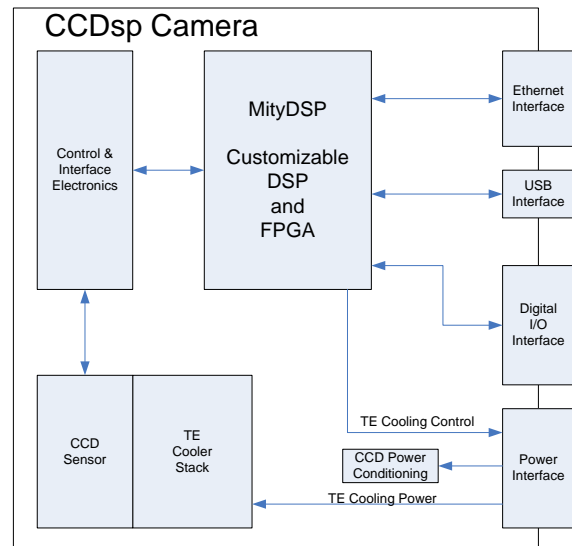


Figure 2: MityCCD Block Diagram

TECHNICAL SPECIFICATIONS

A summary of the imaging performance of the MityCCD-H10141 family of cameras is included below. This family of cameras utilizes sensors from the Hamamatsu S10141 family of CCDs. Detailed datasheets regarding the sensor performance curves for sensitivity, well depth, quantum efficiency and dark current, etc. may be located on the manufacturer's website [1].

PERFORMANCE

	Min	Typical	Max	Units
Active Pixels	-	1024x122	-	H x V
Pixel Size	-	12 x 12	-	W x H ; μm
Imaging Area	-	12.28 x 1.46	-	W x H ; mm
Gain, Setting 1 ²		4.7		e ⁻ per count
Gain, Setting 2 ²		14.2		e ⁻ per count
ADC resolution	-	16	-	bits
Readout Noise @ 500 kHz, Gain 1 / 2	-	12 / 22	18 / 30	e ⁻
Dynamic Range, Gain 1 / 2	-	27K / 44K	-	
Readout Rate ¹	20	493	500	kHz
Full Frame Rate	-	3	-	Frames / s
CCD Binned Frame Rate ⁴	-	200	-	Frames / s
USB Link Throughput ³	-	13,000	-	Kbytes / s
Ethernet Link Throughput ³	-	1,000	-	Kbytes / s
Cooling capability below ambient	35	40	-	°C
Power Utilization ⁵	8	21	31	Watts
Power Input ⁵	12	18	24	V DC

Table 1: MityCCD-H10141 Performance

Note: Data in Table 1 is specific to the MityCCD-H101411007 camera. The Full Frame Rate and CCD Binned Frame Rate would be slower for the MityCCD-H101411008, MityCCD-H101411009, MityCCD-H101411107, MityCCD-H101411108 and MityCCD-H101411109 cameras.

	Typical	Units
Pixel Well Depth, typical	75	k e ⁻ / pixel
Readout Well Depth, typical (for binning)	150	k e ⁻ / pixel
Dark Current @ -15 °C	1.5	e ⁻ / pixel / s

Table 2: S10141 Electrical and Optical Characteristics

Notes:

- 1 – The Readout Rate in a MityCCD stock camera is 493 KHz. The readout rate may be customized for user specific applications, please contact Critical Link, LLC, for details.
- 2 – The MityCCD cameras offset two gain settings, which are user configured via a software setting.
- 3 – Throughput rates assume a back-end PC capable of accepting the data rates provided by the camera.
- 4 – Assumes Full Vertical Binning within the CCD down to 1 output row using USB interface
- 5 – The MityCCD camera must be powered by the MityCCD-PS power supply assembly, which uses a 12-24 DC input in order to generate all necessary voltages to run the camera. The 8 Watt minimum power utilization represents a camera with no TE cooling enabled.

QUANTUM EFFICIENCY

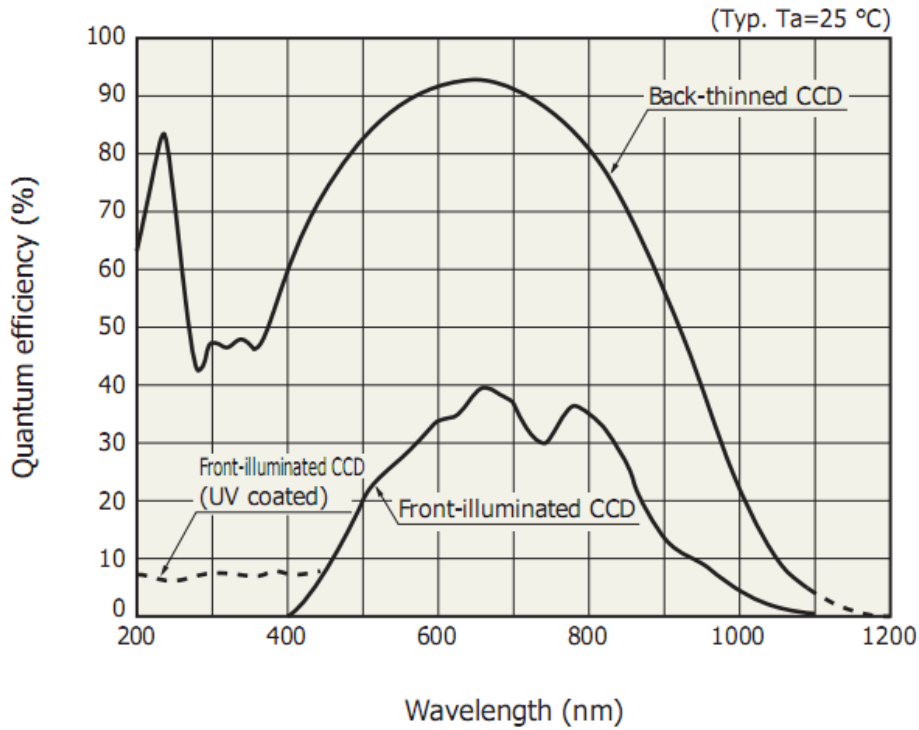



Figure 3: Spectral Response

OPERATING AND STORAGE CONDITIONS

Ambient Temperature Range	0 C to 30 C
Humidity	< 80%, Non-Condensing
Storage Temperature Range	-20 C to 55 C

Table 3: MityCCD-H7031 Operating and Storage Conditions

ORDERING INFORMATION

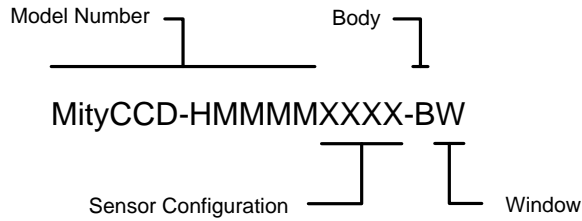

Model Number	Description
MityCCD-H10141	Back-Illuminated

Sensor Config.	Description
1007S	1024 x 122
1008S	1024 x 250
1009S	1024 x 506
1107S	2048 x 122
1108S	2048 x 250
1109S	2048 x 506

Body	Description
A	Standard Body
B	Open Frame – No Flange required
C	Reserved
D	Enclosed Body
E	Reserved

Window	Description
S	AR-coated Sapphire Window

Additional Notes:

1. A MityCCD-H10141 camera also requires a power supply package and flange package that can be ordered separately. For more information contact Critical Link, LLC for details.
2. A MityCCD-H10141 camera also requires a software package. For more information contact Critical Link, LLC for details.
3. A MityCCD-H10141 camera can be ordered with optional accessories. For more information contact Critical Link, LLC for details.

EXTERNAL DIMENSIONS

Unless otherwise noted, all dimensions are in mm [inches].

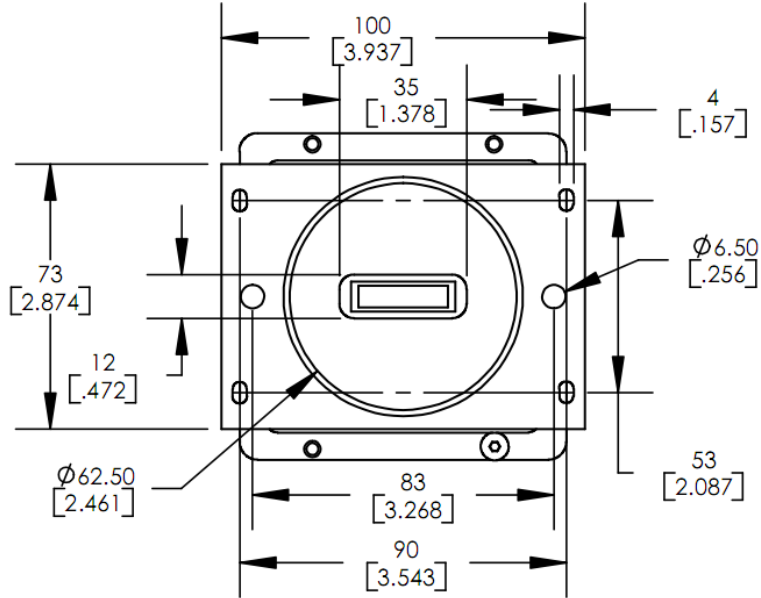


Figure 4: Standard Body (Type A) with Rectangular Flange (Type A) - Front View

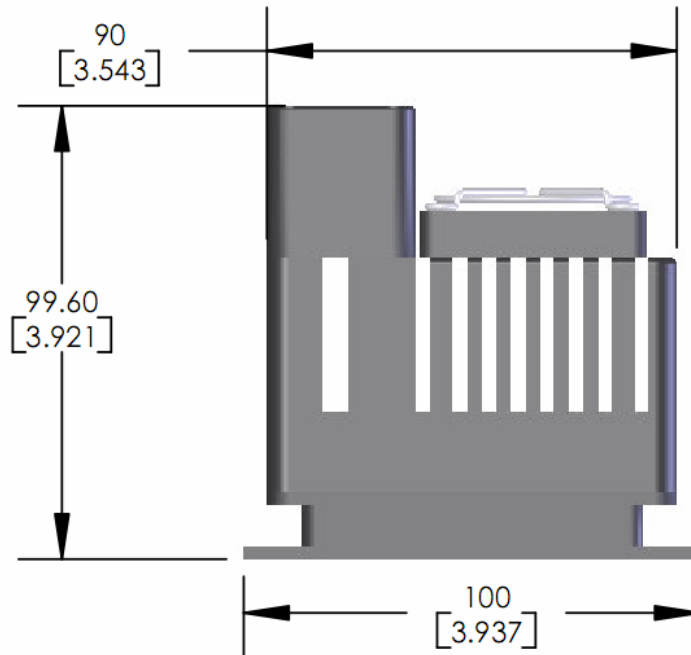


Figure 5: Standard Body (Type A) with Rectangular Flange (Type A) - Side View

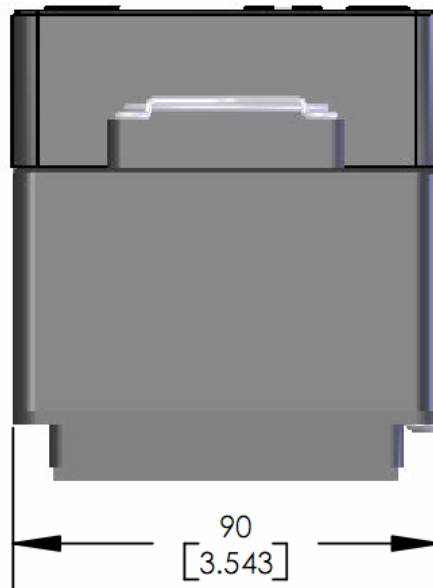


Figure 6: Standard Body (Type A) with Rectangular Flange (Type A) - Top View

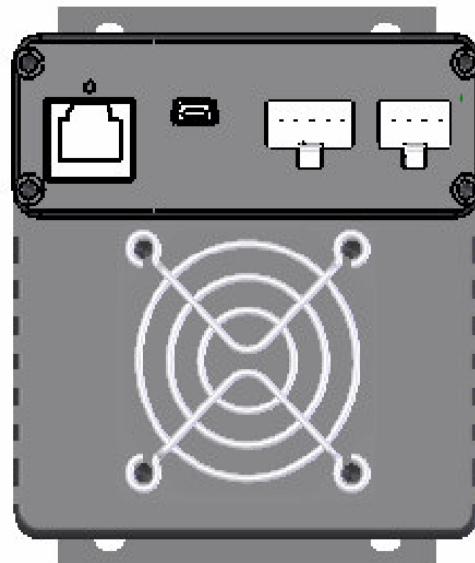


Figure 7: Standard Body (Type A) with Rectangular Flange (Type A) - Back View

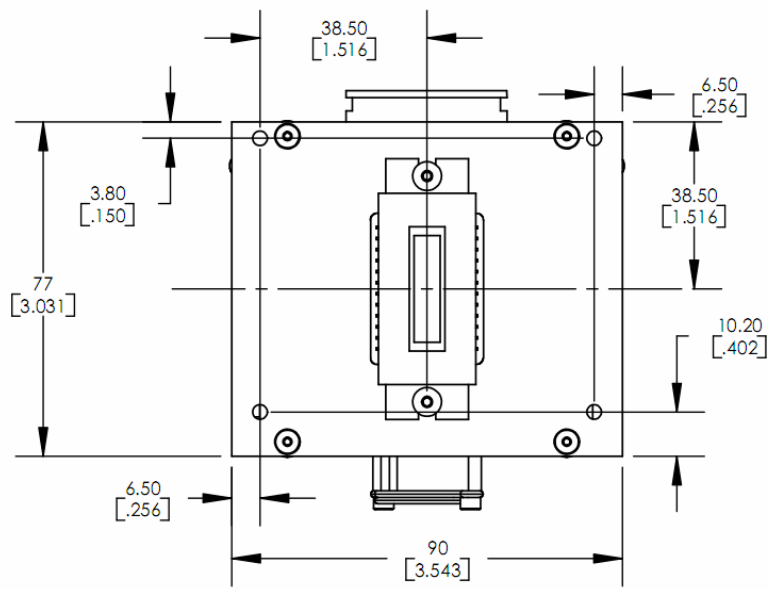


Figure 8: Open Frame Body (Type B) – Front View

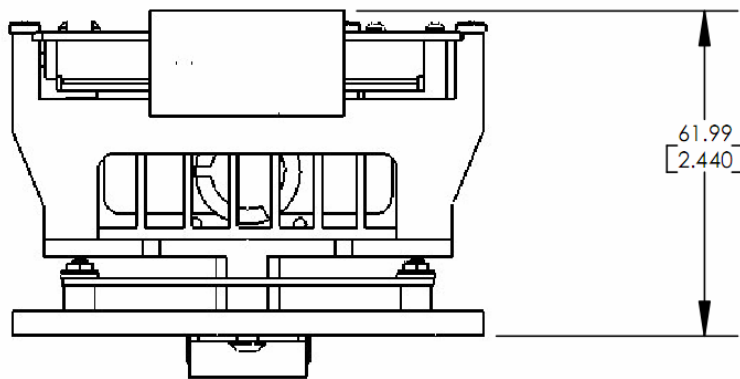


Figure 9: Open Frame Body (Type B) – Top View

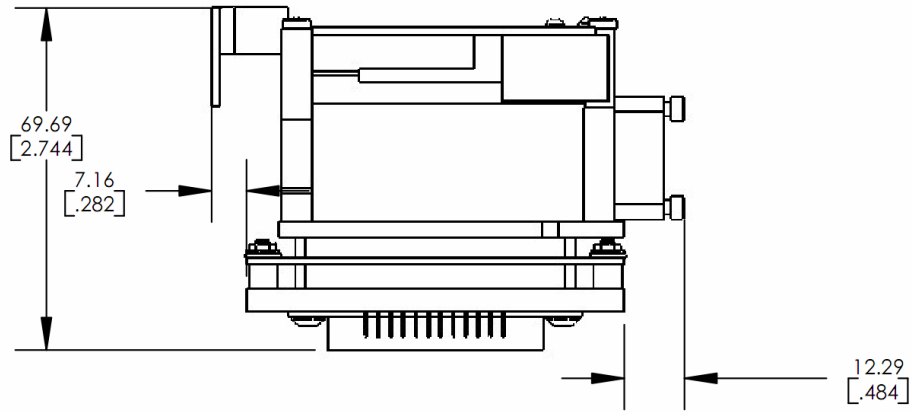


Figure 10: Open Frame Body (Type B) - Side View

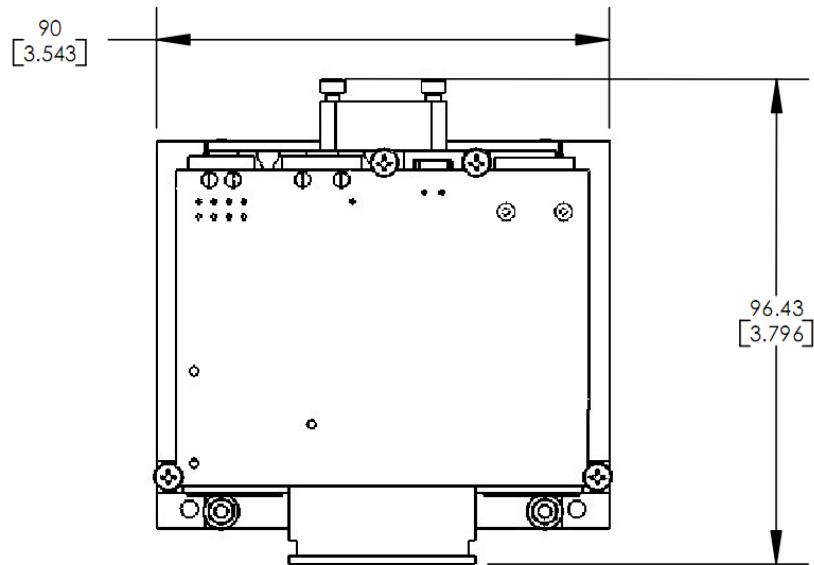


Figure 11: Open Frame Body (Type B) - Back View

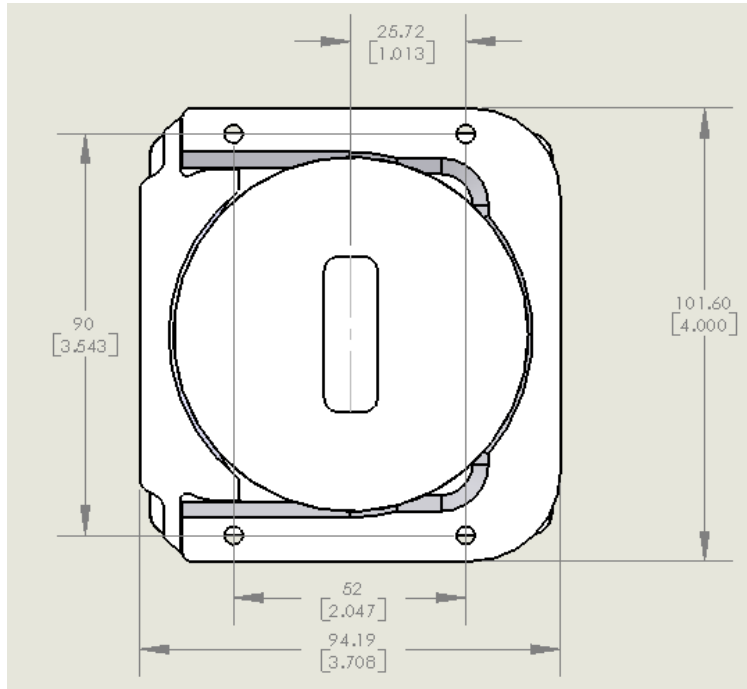


Figure 12: Enclosed Body (Type D) with Split Flange (Type C) – Front View

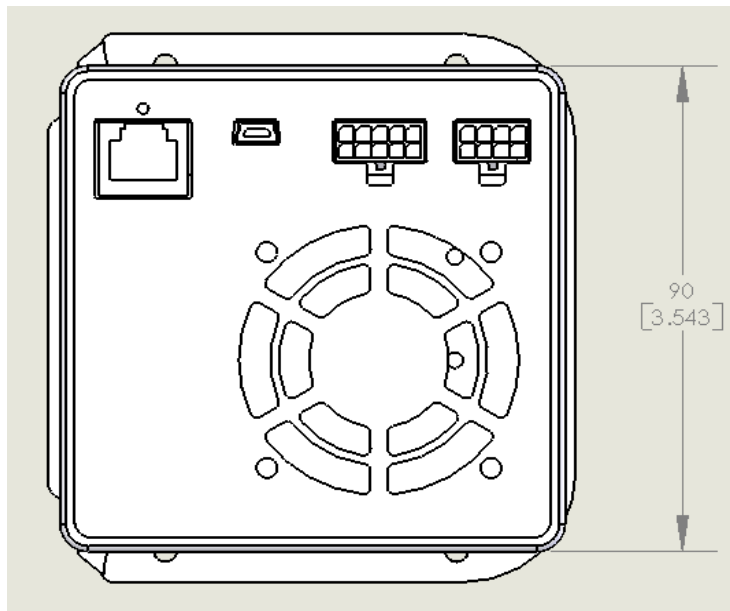


Figure 13: Enclosed Body (Type D) with Split Flange (Type C) - Back View

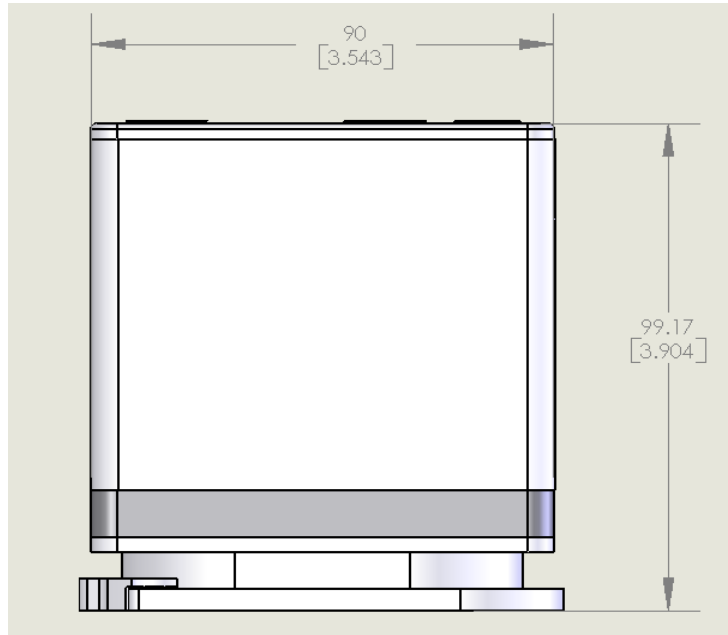


Figure 14: Enclosed Body (Type D) with Split Flange (Type C) - Top View

ELECTRICAL CONNECTIONS

In addition to industry standard Mini USB and 10/100 Mbit based Ethernet RJ-45 communications interface jacks, the MityCCD also provides the following interface connectors.

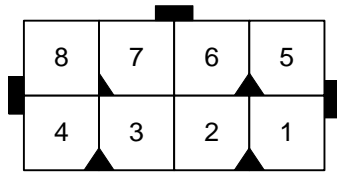


Figure 15: Power Connector¹ - Rear View.

Pin	I/O	Description
1	O	TE Cooler Control
2	I	TE Cooler -V Rtn
3	-	GND
4	I	+6 V (1 Amp Max)
5	I	TE Cooler +V In (7 V at 2 Amps Max)
6	I	+ 15 V (0.3 Amp Max)
7	I	+32 V (0.04 Amp Max)
8	I	-9 V (0.2 Amp Max)

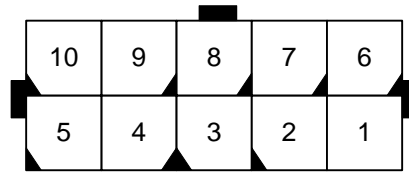


Figure 16: GPIO Connector² - Rear View.

Pin	I/O	Description
1	I/O	General Purpose TTL I/O 3
2	I/O	General Purpose TTL I/O 1
3	O	+5 VD (800 mA max)
4	I	RS232 RX
5	-	GND
6	-	GND
7	I/O	General Purpose TTL I/O 2
8	I/O	General Purpose TTL I/O 0
9	-	GND
10	O	RS232 TX

Notes:

1 – Mating connector is Mfg part # 43025- 0800, please contact Critical Link, LLC, for details.

2 – Mating connector is Mfg part # 43025- 1000, please contact Critical Link, LLC, for details.

REVISION HISTORY

Table 9: Revision History

Date	Rev	Change Description
20-Sep-19	A	Configuration Management release / Corrected read-out rate in table 2

REFERENCE

- [1] Hamamatsu Photonics K.K., Solid State Division, "CCD area image sensor S10140/S10141 series," Jul 2009
http://jp.hamamatsu.com/resources/products/ssd/pdf/s10140-1007_etc_kmpd1094e06.pdf